

## 635: GRACILINEMA MEDIUM

D-Glucose	2.00	g
Tryptone (BD BACTO)	2.00	g
Sodium resazurin (0.1% w/v)	0.50	ml
L-Cysteine HCl x H <sub>2</sub> O	0.25	g
<b>Trace element solution SL-10</b>	10.00	ml
<b>Modified Wolin's mineral solution</b>	25.00	ml
<b>Wolin's vitamin solution (10x)</b>	1.00	ml
Distilled water	970.00	ml

Dissolve ingredients (except cysteine, trace elements and vitamins), sparge medium with 100% N<sub>2</sub> gas for 30 - 45 min to make it anoxic. Then add cysteine, adjust pH to 7.0, dispense medium under same gas atmosphere in anoxic Hungate-type tubes or serum vials and autoclave. After sterilization add trace elements and vitamins from sterile anoxic stock solutions prepared under 100% N<sub>2</sub> gas. Vitamins should be sterilized by filtration. Check pH of complete medium and adjust to 7.0, if necessary.

### Trace element solution SL-10 (from medium 320)

HCl (25%)	10.00	ml
FeCl <sub>2</sub> x 4 H <sub>2</sub> O	1.50	g
ZnCl <sub>2</sub>	70.00	mg
MnCl <sub>2</sub> x 4 H <sub>2</sub> O	100.00	mg
H <sub>3</sub> BO <sub>3</sub>	6.00	mg
CoCl <sub>2</sub> x 6 H <sub>2</sub> O	190.00	mg
CuCl <sub>2</sub> x 2 H <sub>2</sub> O	2.00	mg
NiCl <sub>2</sub> x 6 H <sub>2</sub> O	24.00	mg
Na <sub>2</sub> MoO <sub>4</sub> x 2 H <sub>2</sub> O	36.00	mg
Distilled water	990.00	ml

First dissolve FeCl<sub>2</sub> in the HCl, then dilute in water, add and dissolve the other salts. Finally make up to 1000.00 ml.

### Modified Wolin's mineral solution (from medium 141)

Nitrolotriacetic acid	1.50	g
MgSO <sub>4</sub> x 7 H <sub>2</sub> O	3.00	g
MnSO <sub>4</sub> x H <sub>2</sub> O	0.50	g
NaCl	1.00	g
FeSO <sub>4</sub> x 7 H <sub>2</sub> O	0.10	g
CoSO <sub>4</sub> x 7 H <sub>2</sub> O	0.18	g
CaCl <sub>2</sub> x 2 H <sub>2</sub> O	0.10	g

ZnSO <sub>4</sub> x 7 H <sub>2</sub> O	0.18	g
CuSO <sub>4</sub> x 5 H <sub>2</sub> O	0.01	g
AlK(SO <sub>4</sub> ) <sub>2</sub> x 12 H <sub>2</sub> O	0.02	g
H <sub>3</sub> BO <sub>3</sub>	0.01	g
Na <sub>2</sub> MoO <sub>4</sub> x 2 H <sub>2</sub> O	0.01	g
NiCl <sub>2</sub> x 6 H <sub>2</sub> O	0.03	g
Na <sub>2</sub> SeO <sub>3</sub> x 5 H <sub>2</sub> O	0.30	mg
Na <sub>2</sub> WO <sub>4</sub> x 2 H <sub>2</sub> O	0.40	mg
Distilled water	1000.00	ml

First dissolve nitrilotriacetic acid and adjust pH to 6.5 with KOH, then add minerals. Adjust final to pH 7.0 with KOH.

### Wolin's vitamin solution (10x) (from medium 120)

Biotin	20.00	mg
Folic acid	20.00	mg
Pyridoxine hydrochloride	100.00	mg
Thiamine HCl	50.00	mg
Riboflavin	50.00	mg
Nicotinic acid	50.00	mg
Calcium D-(+)-pantothenate	50.00	mg
Vitamin B <sub>12</sub>	1.00	mg
p-Aminobenzoic acid	50.00	mg
(DL)-alpha-Lipoic acid	50.00	mg
Distilled water	1000.00	ml