

718a: PETROTOGA HALOPHILA MEDIUM

KCl	0.34	g
MgCl ₂ x 6 H ₂ O	4.00	g
MgSO ₄ x 7 H ₂ O	3.45	g
NH ₄ Cl	0.25	g
CaCl ₂ x 2 H ₂ O	0.14	g
K ₂ HPO ₄	0.14	g
NaCl	50.00	g
Modified Wolin's mineral solution	10.00	ml
Fe(NH ₄) ₂ (SO ₄) ₂ x 7 H ₂ O (0.1% w/v)	2.00	ml
Sodium resazurin (0.1% w/v)	0.50	ml
NaHCO ₃	1.00	g
Yeast extract	2.00	g
D-Glucose	5.00	g
Wolin's vitamin solution (10x)	1.00	ml
Na ₂ S x 9 H ₂ O	0.50	g
Distilled water	1000.00	ml

Dissolve ingredients (except bicarbonate, yeast extract, glucose, vitamins and sulfide), then sparge medium with 80% N₂ and 20% CO₂ gas mixture for 30 - 45 min to make it anoxic. Add bicarbonate, dispense medium under same gas atmosphere into anoxic Hungate-type tubes or serum vials and autoclave. After autoclaving add yeast extract, glucose, vitamins and sulfide from sterile anoxic stock solutions prepared under 100% N₂ gas. Vitamins should be sterilized by filtration. Prior to inoculation adjust pH of complete medium to 6.5 - 6.7, if necessary.

Modified Wolin's mineral solution (from medium 141)

Nitrilotriacetic acid	1.50	g
MgSO ₄ x 7 H ₂ O	3.00	g
MnSO ₄ x H ₂ O	0.50	g
NaCl	1.00	g
FeSO ₄ x 7 H ₂ O	0.10	g
CoSO ₄ x 7 H ₂ O	0.18	g
CaCl ₂ x 2 H ₂ O	0.10	g
ZnSO ₄ x 7 H ₂ O	0.18	g
CuSO ₄ x 5 H ₂ O	0.01	g
AlK(SO ₄) ₂ x 12 H ₂ O	0.02	g
H ₃ BO ₃	0.01	g
Na ₂ MoO ₄ x 2 H ₂ O	0.01	g
NiCl ₂ x 6 H ₂ O	0.03	g
Na ₂ SeO ₃ x 5 H ₂ O	0.30	mg
Na ₂ WO ₄ x 2 H ₂ O	0.40	mg

Distilled water	1000.00	ml
-----------------	---------	----

First dissolve nitrilotriacetic acid and adjust pH to 6.5 with KOH, then add minerals. Adjust final to pH 7.0 with KOH.

Wolin's vitamin solution (10x) (from medium 120)

Biotin	20.00	mg
Folic acid	20.00	mg
Pyridoxine hydrochloride	100.00	mg
Thiamine HCl	50.00	mg
Riboflavin	50.00	mg
Nicotinic acid	50.00	mg
Calcium D-(+)-pantothenate	50.00	mg
Vitamin B ₁₂	1.00	mg
p-Aminobenzoic acid	50.00	mg
(DL)-alpha-Lipoic acid	50.00	mg
Distilled water	1000.00	ml