

**1167. NSY-MEDIUM FOR POLYNUCLEOBACTER**

Nutrient broth	1.0	g
Peptone from soybean (e.g. Bacto Soytone)	1.0	g
Yeast extract	1.0	g
Tap water	1000.0	ml

The following inorganic solution may substitute for tap water:

MgSO <sub>4</sub> x 7 H <sub>2</sub> O	75.0	mg
Ca(NO <sub>3</sub> ) <sub>2</sub> x 4 H <sub>2</sub> O	43.0	mg
NaHCO <sub>3</sub>	16.0	mg
KCl	5.0	mg
K <sub>2</sub> HPO <sub>4</sub> x 3 H <sub>2</sub> O	3.7	mg
Trace element solution SL-4 (see DSM-medium 14)	0.1	ml
Distilled water	1000.0	ml

Adjust to pH 7.2.