

894. BENNETT'S MODIFIED MEDIUM (P. AGRAWAL, UNPUBLISHED)

Lab lemco	10.0	g
Yeast extract	2.0	g
Glucose	10.0	g
Starch	100.0	mg
Tryptose	2.0	g
CaCO ₃	100.0	mg
CoCl ₂	traces	-
Ferric ammonium	traces	-
Agar	15.0	g
Distilled water	1000.0	ml

Adjust pH to 7.0 and sterilise at 121°C for 15 minutes.